



Clean Water. Fertile Soil. Serious Science.

334 Hecla Street, Lake Linden, MI 49945 USA

Toll Free: 1.888.NITRATE
Orders: 1.906.296.1115
Tech Support: 1.906.296.1130
Fax: 1.906.296.8003

info@nitrate.com
sales@nitrate.com
tech@nitrate.com
www.nitrate.com

Nitrate Reductase Reagent Packs for Automated Laboratory Equipment

Compatible with High Capacity Discrete Analyzers (Thermo Scientific (Konelab) Arena and Aquakem, etc.)

SKU: *NRPk-HDA-3At* (Previously *DA-ARK-1*)

Introduction

Nitrate Reductase Reagent Packs for Discrete Analyzers were developed by NECi in conjunction with the USGS to replace toxic cadmium in automated colorimetric nitrate determination methods. All reagents for the nitrate reductase method are water soluble and stable in solution, making this method optimal for use in discrete analyzers². NECi's Nitrate Reductase from *Arabidopsis thaliana* (AtNaR) is a recombinant enzyme produced by strict QA/QC procedures with reagent grade water, sugars, and salts to ensure lot-to-lot consistency, reliability, and stability. Nitrate Reductase is non-toxic and environmentally benign.

Method

Nitrate Reductase catalyzes the reduction of nitrate to nitrite with the natural electron donor, NADH (reduced nicotinamide dinucleotide) to drive the conversion¹. Nitrate is then reacted with sulfanilamide to form a diazo compound which is then reacted with N-(1-naphthyl)-ethylenediamine dichloride to produce a pink color. This color is then quantitatively measured using a colorimeter or spectrophotometer reading absorbance at 540 nm \pm 20 nm. Results are then analyzed by generating a standard curve using a spreadsheet program or statistics package.

Reagents Included

This Reagent Kit contains **Nitrate reductase from *Arabidopsis thaliana* (AtNaR)**, enzyme diluent, NADH, and instructions:

- AtNaR – 3 units/vial (freeze-dried) in vacuum sealed foil pouch
- Enzyme diluent – 1 squeeze bulb (pre-measured) with reconstitution instructions
- NADH – 2 mg (dry powder in vial) in vacuum sealed foil pouch

Preparation of Additional Reagents and Protocol for the Thermo Scientific Aquakem are explained in detail on the reverse of this sheet

Instrument

This kit (*NRPk-HDA-3At*) is designed for the Thermo Scientific (Previously Konelab) Aquakem³ and Arena, and is also compatible with other high capacity discrete analyzers with similar specifications. Reagent Packs with detailed instructions are also available for other discrete analyzers, flow injection analyzers, continuous and segmented flow analyzers, and other automated laboratory instruments.

Please inquire about adapting our method for other automated laboratory instruments or ask your sales representative.

References

1. Campbell, Wilbur H., GG Barbier, P Song (2006a) Nitrate Reductase for Nitrate Analysis in Water. Environmental Chemistry Letters 4:69-74.
2. Campbell, Wilbur H., Ellen R. Campbell, Lynn Egan (2006b) Green Chemistry Nitrate Determination: An Alternative Nitrate Analysis Method. American Laboratory, February, 2006.
3. Campbell, Wilbur H., ER Campbell (2007) Nitrate Analysis using Different Nitrate Reductase Isozymes. American Lab Aug07 pp45-46.
4. Patton, C. J., et al. (2008) Nitrate Analysis with Nitrate Reductase on the Discrete Analyzer.

Preparation of Buffers and Reagents (not provided in pack):

- **Ethylenediamine tetraacetic acid (EDTA, 25 mM):** Dissolve 9.3 g Ultrapure EDTA (FW = 372.24) in approximately 800 mL deionized water (DI water) contained in a 1 L volumetric flask. Dilute to the mark with DI water and mix well. Transfer to a bottle and store at room temperature. Stable for one year.
- **Phosphate Buffer (pH = 7.5):** Dissolve 3.75 g potassium di-hydrogen phosphate (KH₂PO₄, FW = 136.1) and 1.4 g potassium hydroxide (KOH, FW = 56.11) in about 800 mL of DI water contained in a 1 L volumetric flask. Add 1 mL of the 25 mM EDTA and dilute the resulting solution to the mark with DI water and mix it well. Transfer this solution to a bottle where it is stable at room temperature for about 1 year.
- **Sulfanilamide (SAN):** Pour 300 mL of concentrated HCl into about 500 mL of deionized water contained in a 1 L volumetric flask. Swirl the flask to mix its contents. Dissolve 10 g SAN (FW = 172.2) in the resulting HCl solution, dilute it to the mark with DI water, and mix it well. Transfer this solution to a bottle where it is stable at room temperature for at least six months.
- **N-(1-Naphthyl)ethylenediamine dihydrochloride (NED):** Dissolve 1 g of NED (FW = 259.2) to about 800 mL of deionized water contained in a 1 L volumetric flask. Dilute the resulting solution to the mark with DI water and mix it well. Transfer this solution to a brown glass bottle where it is stable at room temperature for at least six months.

Preparation of Enzyme Reagents (AtNaR and NADH):

1. Reconstitute AtNaR enzyme [clear vial, red cap] using Enzyme Diluent supplied with the kit (detailed instructions included in pack). This gives you a Stock Solution of 3.0 Units/ml. Store this solution in a freezer between uses. Enzyme diluent contains glycerol and other protein stabilizers. **Never freeze enzyme solution in buffer alone! Use the diluent!**
2. To make sufficient reagents for 8 hours of run time, or ~330 assays, transfer contents of the AtNaR Stock Solution to 19 ml Phosphate Buffer (as prepared above) for a total volume of 20 ml.
 - a. For fewer samples or shorter run times, use smaller volumes (e.g. 335 µl AtNaR Stock Solution and 6.3 ml Phosphate buffer for 100 assays in 2.5 hr).
3. Add 1.0 ml Phosphate Buffer to NADH vial [2.0 mg NADH]. For 8 hr/330 assays, dilute to 10 ml total volume using Phosphate Buffer (above)
 - a. For 100 assays/2.5 hr, add 335 µl of this NADH solution to 3.0 ml buffer. Store unused reagent in a freezer.

Protocol for preparing your own NADH and Nitrate Reductase Reagents:

- **NADH Stock (Sigma #N-8129):** Dissolve 0.100 g of NADH (FW = 709.4) in approximately 40 mL of DI water contained in a 50-mL volumetric flask (final concentration 2 mg NADH/ml). Dilute to the mark and mix well. Transfer 1-mL aliquots to 1.5-mL snap-top vials and store at -20°C. Stocks stable for 1 month.
 - a. Working solution: 1 (1 mL) aliquot/9 mL phosphate buffer (10.0 mL total volume).
- **Nitrate reductase from *Arabidopsis thaliana* (AtNaR):** Add 1 mL of phosphate buffer to a 3.0-unit vial of AtNaR. Gently invert the vial several times over the course of at least 10 minutes to affect dissolution. Add the reconstituted enzyme to 19 mL phosphate buffer (20 mL total volume). Rinse the vial several times with aliquots of the final enzyme/phosphate buffer solution. After freeze-dried AtNaR is reconstituted in phosphate buffer, it is stable for about 8 hours at 4°C. This solution is sufficient for about 330 analyses.

Reaction Conditions for the KoneLab AquaKem/Arena:

1. 55 µL AtNaR Solution
2. 5.0 µL Sample. Mix.
3. 15 µL NADH. Mix.
4. Incubate 600 s at 37°C
5. Add 25 µL SAN. Mix and incubate 120 s at 37°C.
6. Add 25 µL NED. Mix and incubate 120 s at 37°C.
7. Measure absorbance at 540 nm, using secondary wavelength correction at 700 nm.

To adapt this assay to other instruments, modify volumes as required, keeping the above ratios equivalent. Contact NECi directly for assistance modifying this method to other automated laboratory instruments or to build a custom reagent pack at no extra cost.